

Horizon Software



Materials testing, analysis and reporting software






Horizon materials testing software

Tinius Olsen is proud to introduce the next evolution of testing software with our Horizon package. As part of our development process, we have taken the best features of our existing software offerings, including Test Navigator, QMat, and EP600 software, added a host of report writing and data manipulation capabilities and, in the process, created a new, unparalleled testing platform that will make easy work of your materials testing programs, whether they're designed for the demanding rigors of R&D or the charting and analysis functions of QC testing.

One of the first features you see within the Horizon software is its use of the most current Windows environments. These familiar formats make it easy to use and learn, especially since the same familiar functionality is maintained throughout the program.

KEY FEATURES

- Test method library
- Test editor
- Tabbed test and recall area
- Multiple machine control
- Output editor
- Multilingual
- Method editor
- Result editor
- Multifaceted security
- Touchscreen enabled*
- Webcam functionality**
- Report consolidation
- Data importing
- * Requires Windows 8 and a touchscreen.
- ** Requires webcam and support arm

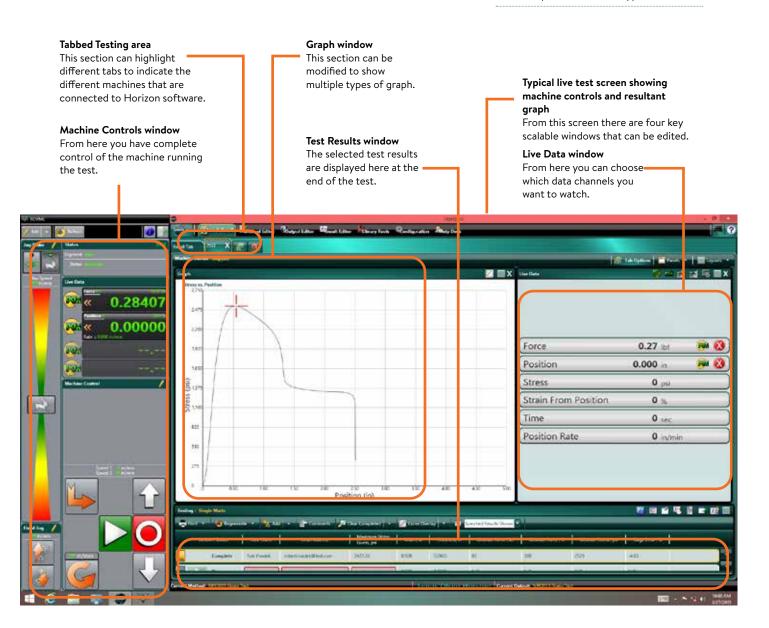




Method library and test area

If your testing regimen follows a quality control analysis to a variety of international standards, then the Test Method Library is probably the first place you will want to visit. Here you can select your desired test methoc which has been written in accordance with different international test standards.

For example, if you need to test for the mel flow index of incoming resin to ASTM D1238 procedure B, you only need to select that routine from the list in the library. Similarly, if you need to determine the pipe crush strengt in accordance with EN802, the tensile strengt of a steel reinforcement bar in accordance with ASTM E8, or the Vicat softening temperature of plastics in accordance with ISO 306, simply make the selection from the library, confirm machine and specimen parameters and start the tests – just four clic of your mouse from start-up to testing!

If, on the other hand, you want to develop your own calculated result from a test, using a national standard as a template, you only have to add the result (and calculation) to the output and save the test set-up with your unique name.

Alternatively, you can develop your own test method to give you complete control over how the test machine performs over the course of the test. You can program the control segments, control options, specimen parameters, report output and how the test machine and software communicate with each other.

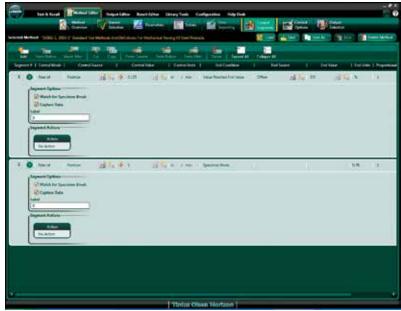

	🔏 🛶 👰 them 🛵 tak									
arent Seath Note ++										
• Open for Switch Query •										
Options	Mitted	L hate	1 Mittodili I							
U Shaw Overviewe	Composition - Familya Position	Entryie Comprised		Compression						
Arthod Hiter	COV 8N 18902 1	CEN (N 1009-1	- A -	Tanvie						
est Type Al	00004 Tear Resistance (Graves Tear) of Plantic First and Sheeting	01004	- 28	Tansle						
article found	DU218 - Met John Text Projecture A	21316	10	Merbuleus Rr						
Bethods Foard	DLDB - Wet lider Test Protecture 8	T1298	112	Wet Indexe Pro						
Otherwy of Working Methods	D1228 Fair Dropepal on Neostance	CORNE	142	Terste -						
	D412 Plantca Texaile - Straw From Ponterr	DATE	23	Terste						
a fartting	DSM Parties Teache - Marin Form Laborationeter	0634	26	Texale						
The base beatly	DER Parties Tessile - Stan Free Poster	D638	14	Taribe						
And instant from the	DIS75 Partici Comprission	10095	15	Compression						
Contraction of the local division of the loc	D Y/D Resure - Shain From Postson	0790	35	Fewers						
Crossel Free III/IIa	0002 Tample Properties of This Plastic Sheving	2882	22	Tansle						
Contractor 1	38 Metals Terrale - 02% Offset, Stain From Externation	100	1.98	Tangle						
and the later of t	Hi Metah Tensle - 52% CPS, 52% Rul, Train Pose Entersconeter	18	14	Tarraile						
The second second	All Merail Tanale - Horsenal VTM	- 11	-10	Tansile						
Contraction of the local division of the loc	164540 Metals Tensle - 82% OF5 IEEE U.S. + Value Street New Extensionater	10154	41	Tamala						
CONTRACTOR DESCRIPTION	13 Metals Compression	18	42	Corposini						
Contraction of the local division of the loc	Bit NC 13854-11899 Meximum Free & Rengelian - Brig Method	#N185138345	45	Territe .						
and the second se	BLIDD 23656-2:389 Maximum Porce - Grab Method	8N 500 18899-2	1.44	Terote						
Transfer to Library	SID 1131 - Met Index Test	80 113	10	Mart Induser Tes						
Characy of Standard Methods	000 527 Plentos Terrale - Share From Extensioneter	80321	-10	Terate						
S carsy or surged without	100-107 Factura Terrate - Strait From Position	80.571	47	Territe						
Tende Sheter	Tatala - Fora va. Reation	Europa Tanalia	18	Tanaba						

Library search result Looking for a standardised tensile test routine.

KEY FEATURES

- Searchable database of international test standards.
- Standards from ASTM, ISO, EN, BS DIN and many more.
- Tests for tension, flexure, melt index, compression etc.
- Ability to customize the test setup using a standard as a template.
- Setups are available in multiple languages and dialects – even create your own!
- Multiple levels and types of security to protect data and equipment.

Control segments setup within the Method Editor section of Horizon software The number of control segments available for each test and test type is unlimited, although experience indicates that typically no more than five are generally used.



Tabbed test and recall area

Test screen showing the tab labeled for multiple MP1200 melt Indexers The software is ready to start the tests on multiple MP1200 or MP600 models etc, while at the same time controlling the tests being performed on the first tab. This example is an H5kT tensile tester.

The next feature you may notice is that Horizon can perform multiple tests at one time, controlling and gathering test data from multiple machines (assuming that your PC has the necessary hardware to control multiple machines). This is true whether you're controlling and gathering data from multiple melt indexers, hydraulic tensile testing machines; and/or electromechanical testing machines - whether they're performing tensile, compression, flexural, tear, peel, or other tests.



KEY FEATURES

- Examine test results from previous tests while performing live tests.
- Ability to test multiple machines and machine types, with sufficient PC power and bandwidth.
- Multiple graph types can be used for each test.
- Full PC networking integration and backup across a PC network.
- Recall function enabling users to add key data perhaps not available, or missed, pre-test.

International State Stat	art Test
134 134 256 266 267 268 268 268 268 268 268 268 268 268 268	
TII Side Tree	100 v. te
	1.00
 In the tas and ta	
Man Temperature 1 cl	100 T
and the second se	- Bank
A Million I and an Annual A	ini. 164
Inter Party Sets 1. Martine N. 1. Annual Inter Database 1.	1 114-
Design 18888 - Design 10	193 BB.2 T
Dates hence prove () park on () 7451 And Dimpt N Indiant set indiant hence Indian hence	
	3 1.07-
T AREAL THE THE THE THE AND THE ADDRESS OF THE ADDR	

Results

A six-station Deflection Temperature Under Load test run on an HDTM603.




Result editor and output editor

Multiple, different graphs can be produced per test, using different measurement axes.


Once all the data has been gathered, Horizon can consolidate it into reports that you can customize to your or your customer's individual needs. The output editor allows unprecedented formatting of your data. You can select what live data can be shown during the test; the acceptable limits of the results; the graphical representation of the test, in multiple formats; the layout of the report including the use of your or your customer's logo; and export or convert the resultant data into another format

These reports can be distributed across one PC, multiple PCs, or even across a network; the presentation of these test reports is compatible with multiple common formats, including ERP.

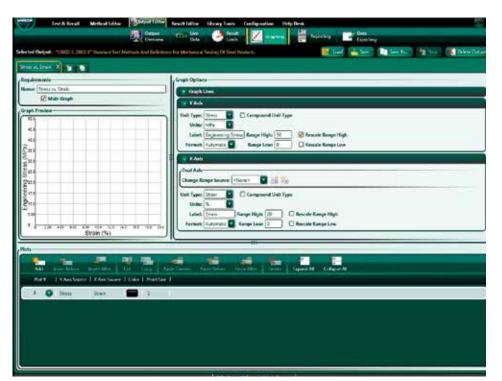

Graph set-up In the Output Editor function of Horizon software.

Customized formulae can be written to create your own unique results In this particular example, the result calculation is a 'standard' result and consequently uneditable and grayed out.

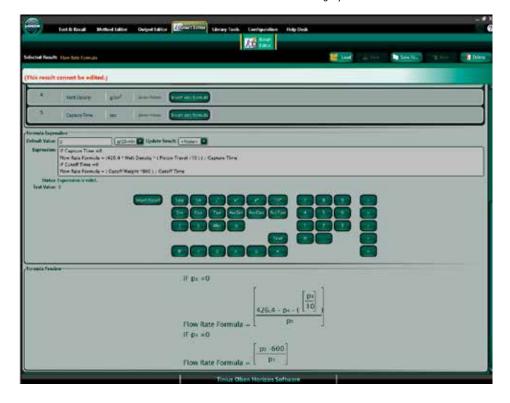

KEY FEATURES

- Multiple graph types can be used per test, such as stress v strain or load v time.
- Unique results can be calculated that draw on parameters and results from other tests and equipment.
- Reports can be generated using your customer's logo.
- Reports can be created in the precise format you desire.
- Data can be exported to a variety of different formats.
- An overview enables you to preview what has been created.



Report generation

	Constra Ba		Law providence		-	_	_	_	
			12.				-		
	Kolona (Janun, Innes Compose 3	in Party in Sec.		Report	Country Seguet	Section limits			Case, Maser, Classow,
Ten ten	E farigen								
🥙 « 0.14	frag bitter								
0.0000	the building has	a de la companya de l	Internet in the		And I have a	Courses and a provider		and in the second second second	
and the rest	10 mm	A1		German-	(internet)		1945		
		(Å.)		444		-	(H))		
	- + • +++	76 Ta ++		Sylvetin		344	18	14 1	
	4 0 mm/m	di la materi		hallow	100	54	141		
	• @ tostion.	and Tay Services		the land	-	-	- 141		
AN AN AN AN		ANT, testimer		441			(H)	19	
	L			-					,
	3 3 7	8 11 0			and the set of the set	com hi			IN CASE A STREET

Result Screen Tabs allow the generation of a single test report for each specimen, and also a summary test report. On each report, individual results from drop-down lists, including operator generated results, can be selected for display.



Insert your or your


customer's logo As with all

sections, the size and location

of this section is user defined.

KEY FEATURES

- Multiple graph types used in the test can be displayed in the report.
- Drop-down list of standard results, and unique results, can be shown in the reports.
- Reports can be generated using your or your customer's logo.
- Single reports can be generated for each completed test; additionally, summary reports can also be generated.
- Common SPC analyses can be preformed and the graphical results printed.

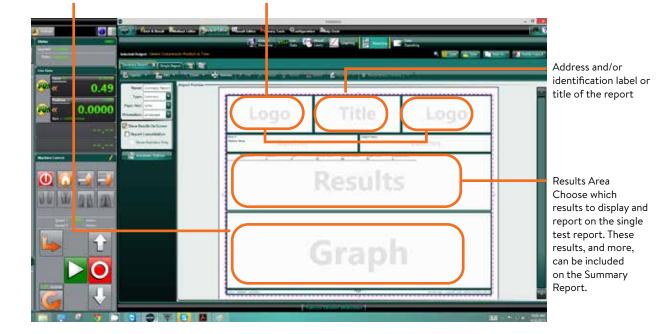

Graph result Set-up allows

axes to be defined along

with the size and location

of the graph(s).



Report generation

HARDWARE REQUIREMENTS

- 2GHz Pentium Dual Core or better.
- 64-bit processor
 - Note: A new installation of Horizon will install with SQL Server 2019 Express by default. SQL Server 2019 Express is NOT supported on 32bit systems.
 - Note: If updating an existing Horizon installation to v10.3.0.0 (or above), a 32-bit system is supported.
- 8GB RAM:
 - Connecting multiple testing machines to a single computer may require additional memory and/or a faster processor.
- 512 MB DirectX 10.0 capable video card or better
- A non-integrated Video card is recommended for best performance.
- 100 GB of available hard disk space (minimum)
 - Note: SQL Server 2019 Express (which installs with Horizon) requires a minimum of 6 GB.
 - Note: Consider more hard drive space for archiving purposes, if testing large quantities of data, or if using a webcam to record videos of tests.
- Machine communication:
 - RS232 Machines (i.e. 398, SR2, S-Series, T-Series, 602, MP993, MP600, Impact and HDV).
 - One port per testing machine is required. The port may be either an integrated RS232 port or USB port using an RS232/USB adapter.
 - USB machines (i.e. MP1200 and VMC controlled).
 - One USB port required per testing machine
- Additional ports:
 - One USB port for use with the software key (required).
 - Ports required for each measuring device, barcode scanner, etc.
 - One USB port and one display port if using touchscreen purchased from Tinius Olsen.
- Mouse and keyboard supported by Windows.
- Monitor: 1600 x 900 (widescreen) or higher.
- Windows-compatible printer, for reporting capabilities.
- Windows-compatible sound card and speakers for audio playback.
- An active internet connection (for TeamViewer use and Help Desk support) is recommended.

SOFTWARE REQUIREMENTS

- A new installation of Horizon v10.3.0.0 (or higher) is designed for a 64-bit Operating System running Windows 10 ONLY.
 - Note: If updating an existing Horizon installation to v10.3.0.0 (or above), windows 7 SP1, 8.1 and 10 are supported.
 - Note: If using Windows 8 (or higher), the '.NET Framework 3.5' Windows feature must be enabled before attempting installation.
- Internet Explorer 8 or higher.
- Once installed, users must be granted full read\ write access to the installed software folders (i.e. the Horizon and Microsoft SQL Server folders).



KEY FEATURES

- Full PC networking integration and backup across a PC network using industry standard SQL database structure.
- Multiple licensing of Horizon for multiple installations to review methods, results and reports away from the testing machine, lab or plant.
- Multi-level user password capability.
- Recall function enables users to add key data perhaps not available, or missed, prior to the test.
- Built in Team Viewer licenses enabling TO engineers to log in directly, upon request, and provide on-machine technical support.
- Live results during test; there is no limit to the number of results displayed and these do not have to be used in the test report.
- With sufficient PC power and bandwidth, multiple testing machines can be run from a single PC.
- Users can use pre-written international standards methods as templates to modify and build their own test methods.



www.tiniusolsen.com info@tiniusolsen.com

Horsham, PA, USA
Redhill, Surrey, UK
Noida, UP, India
Shanghai, PR China